

# **MilkKit**

## **Milk Rapid Test Kit**

## MilkKit Milk Rapid Test Kit

World-class quality at affordable cost



MilkKit tests can be used to monitor antibiotic residues and mycotoxins, melamines, etc in milk and various food samples.

## Current Offer in Dairy Industry

### Veterinary Drug Residue Rapid Test Kit

Major veterinary antibiotics, steroids, hormones, disinfectants, detergent, etc.

### Pesticide Residue Rapid Test Kit

Important pesticides, herbicides, etc.

### Mycotoxin Rapid Test Kit

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
for aflatoxin M1 testing, ranging from EU MRL, USA MRL to Codex MRL, 0.5ppb, 0.05ppb, 0.3ppb, 0.4ppb, etc.

### Milk Nutrition & Milk fraud Rapid Test Kit

for testing of bovine IgG, bovine lactoferrin, vitamin, as well as testing cow milk added into goat or camel milk, etc.


### Test Equipment for Milk Tests

from portable reader system to desktop reader system to ensure the visualize of test data and printing.



# Milk Kit Test
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
## General Step for milk testing



## $\beta$ -lactams & tetracyclines, BT Combo Test Kit

A lateral flow rapid test kit based on receptor assay to detect multiple  $\beta$ -lactams & tetracyclines antibiotics in 10min.

### Ordering information

Product Code	Reaction Mode	Detection Level
TESTMIRA2002A	5'+5', room temp.	EU MRL

.....

## $\beta$ -lactams & tetracyclines & sulfa drugs, BTS TriTest 3in1

A lateral flow rapid test kit based on receptor assay to detect multiple  $\beta$ -lactams & tetracyclines & sulfa antibiotics in 6min.

### Ordering information

Product Code	Reaction Mode	Detection Level
TESTMIRA0015A	3'+3', room temp.	EU MRL

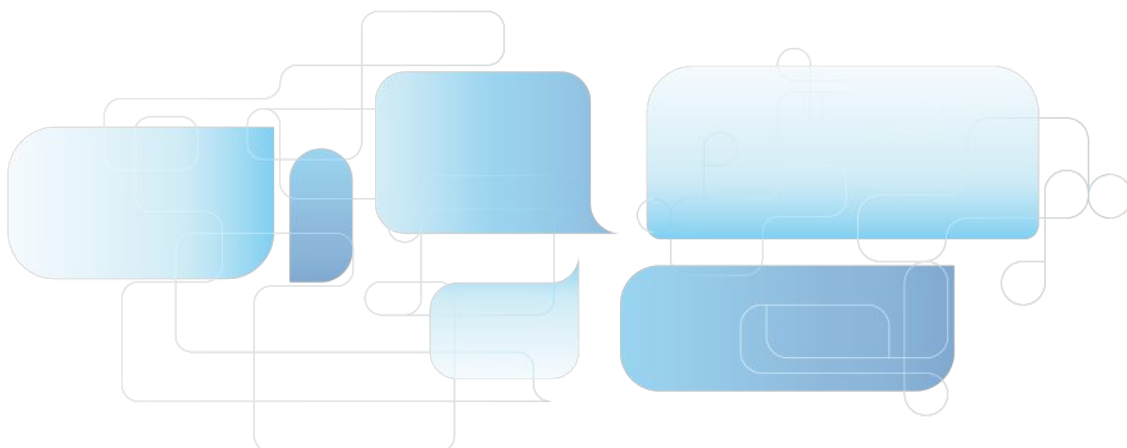
.....

### Kit Component

- A12 tubes of rapid tests, 8 strips per tube
- Microwell holder
- Plastic pipette or micropipette tips
- Kit instruction

### Samples

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.



## β-lactams & tetracyclines & sulfa & quinolones, BTSQ QuaTest 4in1

A lateral flow rapid test kit based on receptor assay to detect multiple β-lactams & tetracyclines & sulfa antibiotics in 6min.

### Ordering information

Product Code	Reaction Mode	Detection Level
TESTMIRA0029A	3'+7', room temp.	EU MRL
.....		

## β-lactams & tetracyclines & streptomycin & CAP, BTSC QuaTest 4in1

A lateral flow rapid test kit based on receptor assay to detect multiple beta-lactams, tetracyclines, streptomycin & chloramphenicol antibiotics in 10min.

### Ordering information

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
Product Code	Reaction Mode	Detection Level
TESTMIRA0018A	3'+7', room temp.	EU MRL
.....		

### Kit Component

- 12 tubes of rapid tests, 8 strips per tube
- Microwell holder
- Plastic pipette or micropipette tips
- Kit instruction

### Samples

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.



## Aminoglycosides, GNKS QuaTest 4in1

A lateral flow rapid test kit based on receptor assay to detect multiple gentamicin, neomycin, kanamycin & streptomycin antibiotics in 10min.

### Ordering information

Product Code	Reaction Mode	Detection Level
TESTMIRA0025A	3'+7', room temp.	EU MRL

.....

## Quinolones, macrolides, lincomycin, erythromycin, QMLE QuaTest 4in1

A lateral flow rapid test kit based on receptor assay to detect multiple quinolones, macrolides, lincomycin, erythromycin antibiotics in 10min.

### Ordering information

Product Code	Reaction Mode	Detection Level
TESTMIRA0024A	3'+7', room temp.	EU MRL

.....

### Kit Component

- 12 tubes of rapid tests, 8 strips per tube
- Microwell holder
- Plastic pipette or micropipette tips
- Kit instruction

### Samples

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.

## Test Equipment and Devices

### Ordering information

Product Code	Product Description
TESTINR1303A	Reader Fast 2 system - the quick and smaller model
TESTINR1304A	Minipipet for test kit, 200ul
TESTINR1305A	Mini Incubator - heating block for rapid tests
TESTINR1307A	ATP fluorescence detector - ATP Reader

### Q3 system Desktop Rapid Test Reader

Ordering information: TESTINR1301A

Weight	3.0kg
Dimension	320mm x 195mm x 156mm
Wavelength	525±30nm
Precision	C.V. ≤3%
Display	6.2Inch LCD
Printer	Built-in thermal printer
Memory	4GB built-in/flash drive



### Fast 2 system Desktop Rapid Test Reader

Weight	1.0kg
Dimension	210mm x 175mm x 75mm
Wavelength	525±30nm
Precision	C.V. ≤2%
Display	3.5Inch LCD
Printer	Built-in thermal printer
Memory	50000 entries

